Flexibler Ausstoßschlauch für Dispenser Varispenser und Varispenser plus/ Flexible discharge tube for dispenser Varispenser and Varispenser plus

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Before initial operation:

Observe safety instructions, exclusions and limitations of use according to the Operating Manuals of the dispenser Varispenser resp. Varispenser plus.

Operating Exclusions:

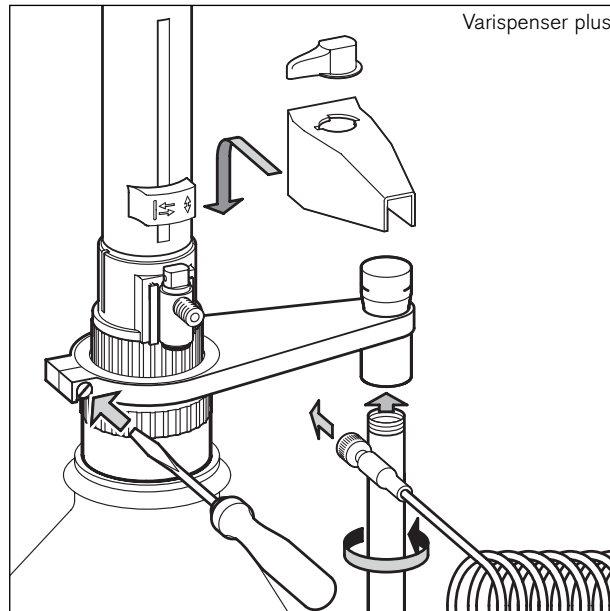
The flexible discharge tube is compatible with all liquids suitable for the dispenser Varispenser or Varispenser plus.

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Montage:

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Attention!

Each time you are going to use the tubing, examine it carefully for any damages. To dispense aggressive liquids, e.g. concentrated acids and lyes, you should take safety measures in addition to the normal precautions of wearing protective clothing, eye protection and protective gloves. We recommend that you use a protective screen.

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
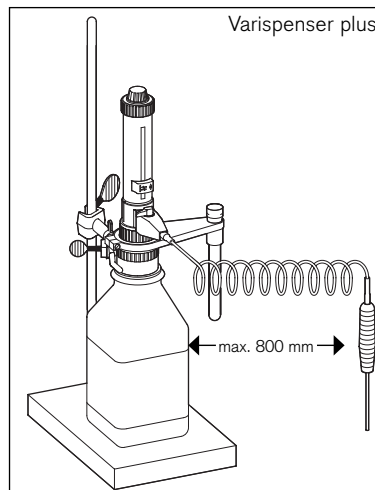
Note:

The coil of the tubing can be stretched to a maximum length of 800 mm. The entire coil must lie in regular loops and must not be twisted.

When working with the flexible discharge tube, the accuracy and coefficient of variation indicated in the Operating Manual are only obtained for volumes ≥ 2 ml and by smooth and steady operation, gently approaching the upper and lower stops. For cleaning rinse the tube carefully. Do not dismantle!

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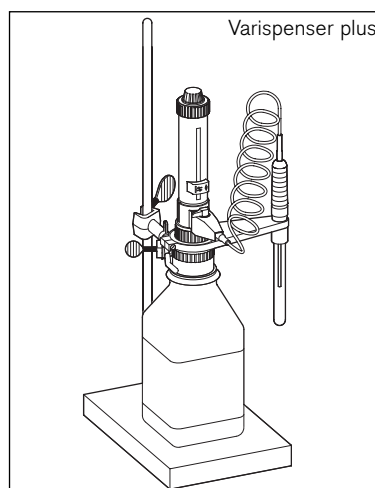Attention!

Use a bottle stand to help prevent the bottle from accidentally tipping over. To help avoid reagent splashing from the tube, always grip the tube firmly by the handle and replace in the holder after use.

Tuyau de d'éjection flexible pour distributeurs Varispenser et Varispenser plus/ Tubo de expulsión flexible para dosificadores Varispenser y Varispenser plus

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