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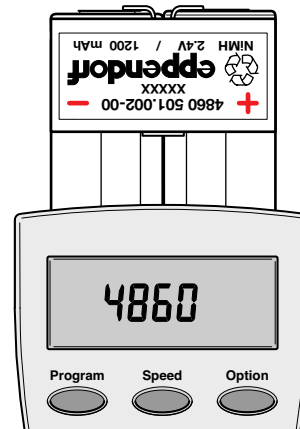
eppendorf® is a registered trademark.
4860 900.101-01/7008 · Printed in Germany

Eppendorf Research® pro

Battery pack
Akku-Pack
Accumulateur
Acumulador
Accumulatore

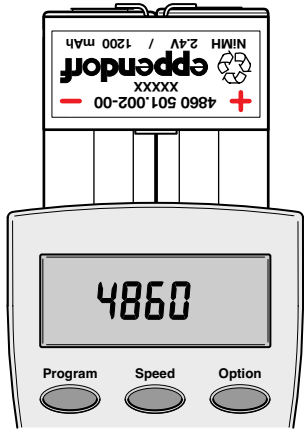
Important information on the Ni-MH 1200 mAh battery

1. Section 3.2 of the instructions describes how to insert the battery.
2. Outside the pipette, the battery poles (+/-) may not be put in contact with metal surfaces.
3. Charge the battery completely before using the pipette for the first time (see Section 3.3 of the instructions). Fully-charged status is indicated by an upright battery symbol in the display as long as the Research pro is located in the charging stand or the charging adapter. During the charging process, the segments of the battery symbol roll continuously.
4. If the Research pro is not going to be used for several months, the battery should be fully charged first and then removed from the battery compartment (observe Item 2!). In order to avoid complete discharge, the removed battery should be recharged approx. every 6 months.
5. Used batteries should be disposed of in accordance with legal regulations.

Batteries should not be disposed of with domestic waste!
Do not heat the battery over 100 °C
and do not throw it on a fire.

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