


Short instructions

Multipette®/Repeater® (X)stream
English (EN)

1 How to use these instructions

These short instructions are a supplement to the operating manual and not a substitute. Read the operating manual before using the Multipette/Repeater (X)stream for the first time.

2 Main illustration

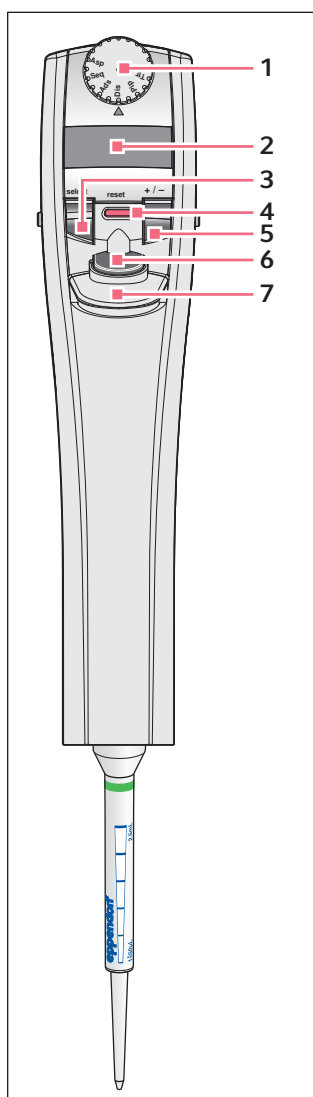

1 Selection dial

Multipette stream/Repeater stream

- *Pip* - Pipetting
- *Dis* - Dispensing
- *Ads* - Auto dispensing

Multipette Xstream/Repeater Xstream

- *Pip* - Pipetting
- *Dis* - Dispensing
- *Ads* - Auto dispensing
- *Asp* - Aspirating
- *Seq* - Sequential dispensing
- *Ttr* - Titration

2 Display

3 Select rocker

For switching between executing and editing parameters.

4 Reset key

For stopping or canceling dispensing operations.

5 +/- rocker

For changing parameters. Higher speed level with volume change when you press the +/- rocker more firmly.

6 Dispensing button

For triggering dispensing operations. For saving entries during editing and for returning to dispensing.


7 Ejector

Releases the inserted Combitip.

Short instructions

Multipette®/Repeater® (X)stream
English (EN)

3 Display

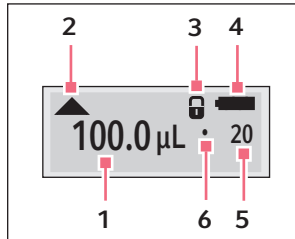

1 Volume

Unit in µl or ml.

2 Direction of the next piston movement.

3 Key lock

🔒 is only displayed if the +/- rocker is locked.

4 Rechargeable battery

Rechargeable battery charging status.

5 Dispensing step

Number of possible dispensing steps.

6 Interpunct

Indicator for the number of dispensing steps.

4 Inserting the Combitip


Keep the ejector pressed down and insert the Combitip straight into the Multipette/Repeater (X)stream.

5 Display after aspirating liquid



After liquid has been aspirated the display shows *Discard next step*. The following step (reverse stroke) must be discarded as the reverse stroke is not a dispensing step.

6 Display after the last dispensing step

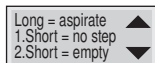

After the last dispensing step you can select between different steps

- *Long = aspirate* - Hold down the dispensing button (> 2 s) - renewed aspiration
- *1 short = no step* - Press the dispensing button once - no function
- *2 short = empty* - Press the dispensing button twice - empty Combitip

7 Ejecting the Combitip

Do not eject the Combitip until it is completely empty.

Your local distributor: www.eppendorf.com/contact

Eppendorf AG · 22331 Hamburg · Germany
eppendorf@eppendorf.com · www.eppendorf.com