

## Installation guide

Software update for BioPhotometer plus 6132

English (EN)

# Software update for BioPhotometer plus 6132

## 1 System requirements

- Windows XP with administrator rights.
- Serial interface (RS232) available on the computer.

## 2 Preparing the devices

Prerequisites


- The BioPhotometer is switched off.
- The computer or the laptop is switched off.

1. Connect the BioPhotometer to the computer or laptop using the serial cable (order no. 6132 869.004).
2. Switch on the computer or laptop.
3. Switch on the BioPhotometer.

## 3 Loading and extracting software

1. Create a folder with a random name on your own computer.
2. Load the file "6131\_V1.37.zip".
  - Download the file from [www.eppendorf-support.com](http://www.eppendorf-support.com).
  - Save the file in the new folder.
3. Extract the file "6131\_V1.37.zip" to the same folder.  
The file "6131\_V1.37.exe" is created in this folder.


## 4 Updating the software

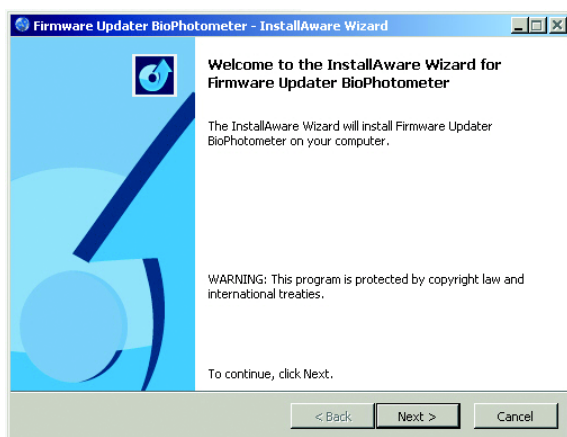


1. Open the file "6131\_V1.37.exe" by double-clicking.  
The Welcome dialog opens.
2. Click *Next* to start the firmware updater.

## Installation guide


Software update for BioPhotometer plus 6132  
English (EN)

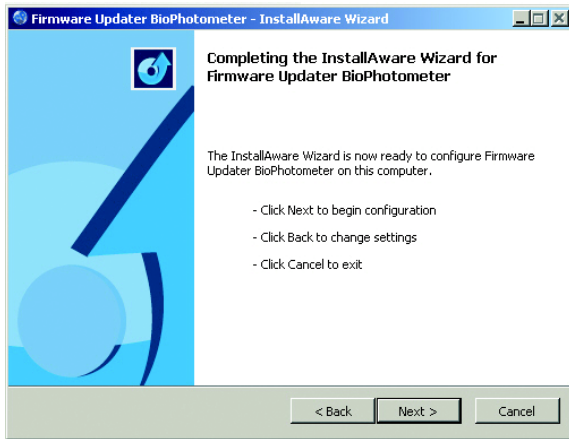

3. Click *Next* to configure the firmware.

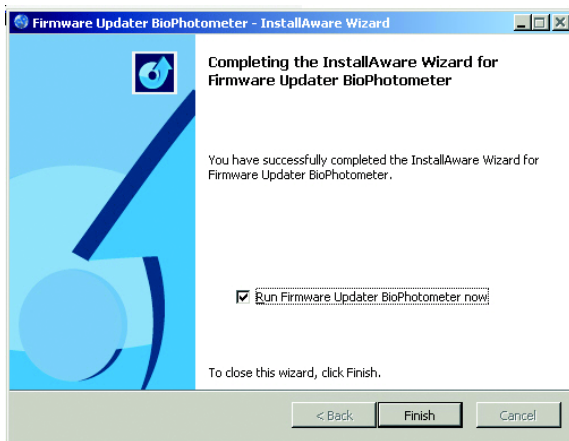

4. Tick the checkbox to start the bootloader.  
5. Click *Finish* to close the firmware updater.  
The bootloader opens.

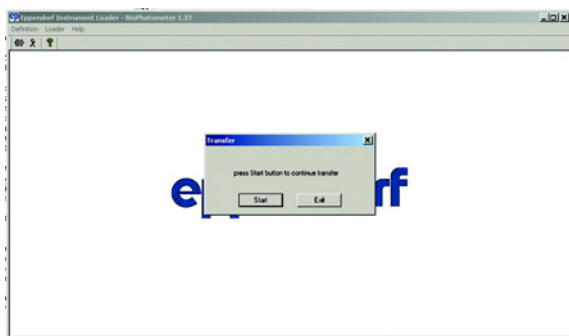

6. Click *Start*.

Your local distributor: [www.eppendorf.com/worldwide](http://www.eppendorf.com/worldwide)  
Eppendorf AG · Hamburg · Germany · Phone: +49 40 538 01-0  
[www.eppendorf.com](http://www.eppendorf.com)