

Update Instruction for Software of Temperature Verification System – Multi-Channel Version 2.1

Requirements

- PC with Windows XP (32 bit)
- MTV Version 1.0 or 2.0 is already installed

Update of application software (for existing systems only)

1. Connect sensor plate to system base (see Operating Manual Temperatur Verification System Multi-Channel).
2. Switch on system base.
3. Connect system base to the notebook (see Operating Manual Temperatur Verification System Multi-Channel).
4. Log in to Windows as Administrator user.
5. Keep MTV installer ready, e.g.
 - MTV.exe on distribution CD
 - MTV.exe on download media
6. Start the MTV installer.



If not stated otherwise, every component shall be installed with its default settings proposed by the installer.

- Ignore potential security warnings through the installation process.
- Make sure that no error message has occurred.

Installation may take some minutes!

7. As soon as installation has finished, start the application software once to actualize the database.

Update Instruction for Software of Temperature Verification System – Multi-Channel Version 2.1

