## Cell culture bottle adapter – Instructions for use

This document supplements the instructions for use of the Centrifuge 5810/5810 R, but does not replace them. Therefore, also read the instructions for use. The current version can be found online at [www.eppendorf.com](http://www.eppendorf.com).

### Product description

The cell culture bottle adapter is designed for the horizontal centrifugation of 25 cm<sup>2</sup> and 75 cm<sup>2</sup> cell culture bottles in the A-4-81 swing-bucket rotor with the MTP/Flex carrier of the Centrifuge 5810/5810 R.

### Safety notes


#### CAUTION! Bottle breakage due to overloading.

- ▶ Never expose the loaded cell culture bottle adapters to more than 1000 x g.
- 
- ▶ Symmetrically load the MTP/Flex carrier of the rotor. When doing so, follow the loading instructions included in the instructions for use of the centrifuge.
  - ▶ Observe the maximum allowed loading height and load (see *Technical data*).
  - ▶ Complete a manual swing-out test when using a bottle type for the first time.

### Centrifugation

1. Place the cell culture bottle adapter in the MTP/Flex carrier so that the seal of the cell culture bottle points inward toward the rotor nut.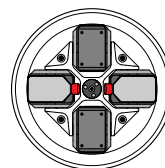

2. Ensure that the cell culture bottle is flush.  
This also applies to the 25 cm<sup>2</sup> cell culture bottle.

Tab. 1: Only the following cell culture bottles\* fit in the cell culture bottle adapter.

75 cm <sup>2</sup> cell culture bottle	Order no.	25 cm <sup>2</sup> cell culture bottle	Order no.
Sarstedt, with filter	83.1811.002	Sarstedt, with filter	83.1810.002
Sarstedt, without filter	83.1811	Sarstedt, without filter	83.1810
		Greiner Bio-One, with filter	690175
		Greiner Bio-One, without filter	690160
		TPP, with filter	90026
		TPP, without filter	90025
		IWAKI, without filter	3102-025

\* ) The centrifugability of the bottles was tested at 32 °C, 1000 x g / 2501 rpm for 90 min.

## Cell culture bottle adapter – Instructions for use

### Cleaning and Disinfection

Use a mild cleaning agent for cleaning.

Use alcohol (ethanol, isopropanol) or disinfectants containing alcohol for the disinfection.

- i** The Cell culture bottle adapter can be autoclaved (121 °C, 20 min).

### Technical data

Max. <i>g</i> -force (rcf)/rotational speed (rpm)	1,000 x <i>g</i> / 2501 rpm *
Max. centrifugation radius	143 mm
Max. load per carrier (adapter, bottle and contents)	380 g
Max. loading height	60 mm
Material	Silicone  > S I <

\*) The cell culture bottles can break at a *g*-force of over 1,000 x *g*. Observe the maximum loading information from the cell culture bottle manufacturer at all times.

### Ordering information

Order No. (International)	Order No. (North America)	Description
5825 719.000	022638972	<b>Adapter</b> for 1 cell culture bottle, used in MTP/Flex buckets, set of 2

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Eppendorf North America, Inc. · 102 Motor Parkway · Hauppauge, N.Y. 11788-5178 · USA