# Exchangeable Thermoblocks for Thermomixers and Thermostats

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**eppendorf**

Copyright© 2012 Eppendorf AG, Hamburg. No part of this publication may be reproduced without the prior permission of the copyright owner.

### **Trademarks**

eppendorf is a registered trademark of Eppendorf AG, Hamburg, Germany.

Thermomixer is a trademark of Eppendorf AG, Hamburg, Germany.

Falcon is a registered trademark of Becton Dickinson, Franklin Lakes, USA.

Registered trademarks are not marked in all cases with <sup>TM</sup> or <sup>®</sup> in this manual

<b>Instructions for use</b> .....	<b>3</b>
<b>Gebrauchsanweisung</b> .....	<b>9</b>
<b>Ordering information/Bestellinformationen</b> .....	<b>15</b>

### 1 User instructions

#### 1.1 Using this manual

These instructions for use for exchangeable thermoblocks are a supplement to the operating manual for the thermomixer. They are not intended as a replacement for this operating manual.

Therefore, also read the operating manual of the corresponding unit before using the exchangeable thermoblock for the first time. The current version of the operating manual can be found in the Internet at [www.eppendorf.com](http://www.eppendorf.com).

### 2 Product description

#### 2.1 Features

The exchangeable thermoblock system for the Thermomixer comfort / Thermomixer R and ThermoStat plus are used for convenient temperature control and mixing of all standard lab tubes. The following exchangeable thermoblocks are available:

Exchangeable thermoblock	Tubes and plates	Use
Exchangeable thermoblock 0.5:	24 × 0.5 ml micro test tubes	<ul style="list-style-type: none"><li>• Mixing and temperature control of the tubes.</li><li>• Use Eppendorf IsoRack to reduce the heat exchange with the environment and maintain the temperature control.</li></ul>
Exchangeable thermoblock 1.5:	24 × 1.5 ml micro test tubes	
Exchangeable thermoblock 2.0:	24 × 2.0 ml micro test tubes	
Exchangeable thermoblock Cryo:	24 × 1.5 or 2.0 ml Cryo tubes	<ul style="list-style-type: none"><li>• Mixing and temperature control of the tubes.</li></ul>
Exchangeable thermoblock lab tubes:	24 tubes at 11 – 11.9 mm, height 30 to 76 mm	

## Instructions for use

Exchangeable thermoblock	Tubes and plates	Use
Exchangeable thermoblock 5.0:	8 × 5.0 ml micro test tubes	<ul style="list-style-type: none"> <li>Mixing and temperature control of the tubes.</li> <li>Mixing possible up to a maximum of 750 rpm.</li> </ul>
Exchangeable thermoblock 15 ml Falcon:	8 × 15 ml Falcon tubes	
Exchangeable thermoblock 50 ml Falcon:	4 × 50 ml Falcon tubes	
Exchangeable thermoblock MTP:	MTP and DWP	<ul style="list-style-type: none"> <li>Mixing and temperature control of the MTP and DWP.</li> </ul> <p>For plates with a membrane bottom, an insertable frame is available which guarantees that the filter does not rest directly on the bottom.</p> <p>DWP cannot always be mixed at the highest frequency (1,400 rpm) as the unit and the plate in the exchangeable thermoblock are stable only at 800 rpm when the total weight is &gt;200 g .</p>
Adapter plate	PCR tubes and plates	<p>An adapter plate is also available for 96 × 0.2 ml PCR tubes or 96-well PCR plates. Max. frequency: 1,300 rpm.</p>
CombiBox	Vials and beakers with heights of up to 60 mm.	<ul style="list-style-type: none"> <li>Only for the ThermoStat plus.</li> <li>Temperature control of vials and beakers up to 95 °C.</li> </ul>

### 3 Intended use

The exchangeable thermoblocks are intended exclusively for indoor use and to hold tubes and plates during sample preparation.

The exchangeable thermoblocks must only be operated by appropriately qualified and instructed personnel. Carefully read the operating manual before working with the exchangeable thermoblocks and familiarize yourself with their use.

### 4 Installation and operation

#### 4.1 Mounting exchangeable thermoblocks

---

**Warning! Injury or property damage from improperly attached exchangeable thermoblocks.**

If the exchangeable thermoblocks are not attached properly, they can become detached from the device or break.

- ▶ Tighten the screws of the exchangeable thermoblocks fingertight.
- 

**Warning! Contamination due to opening tube cover.**

Covers of micro test tubes can snap open involuntarily in the following cases, causing sample material to escape:

- high vapor pressure of the content
  - improperly sealed cover
  - damaged sealing lip
- ▶ Always check that micro test tubes have been sealed tightly before use.
- 

#### 4.1.1 Exchangeable thermoblocks for micro test tubes

Valid for:

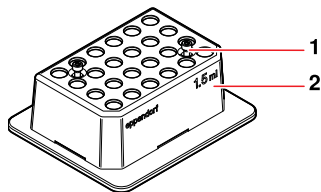
- 0.5 ml micro test tubes
- 1.5 ml micro test tubes
- 2.0 ml micro test tubes
- 5.0 ml micro test tubes
- 11-11.9 mm Diameter of tubes
- 15 ml Falcon tubes
- 50 ml Falcon tubes

### **Warning! Risk of burns from hot surfaces.**

The exchangeable thermoblock can be very hot after heating and cause burns.

- ▶ Allow the heated exchangeable thermoblock to cool down completely before removing the exchangeable thermoblock.

1. Place the exchangeable thermoblock (2) on the thermosocket with the flat edge and the writing facing the front.
2. Tighten the captive screws (1).


### **4.1.2 MTP exchangeable thermoblock**
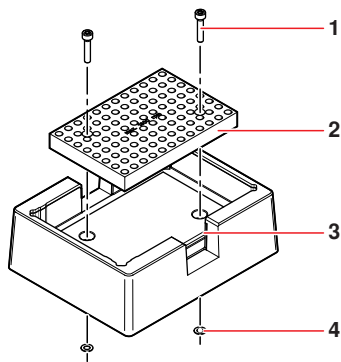
#### **Caution! Melting of the MTP from excessive temperatures.**

- ▶ Only heat the polystyrene MTP up to max. 70 °C.

1. Insert the short screws in the MTP exchangeable thermoblock with the flat washers.
2. Secure the screws from below using the lock washers.  
The lock washers prevent the screws from falling out of the bores if the exchangeable thermoblock is not fixed on the unit.
3. Place the MTP thermoblock on the device in such a way that the spring plate is to the front left.
4. Tighten the screws.

### 4.1.3 Adapter plate for 0.2 ml tubes or 96 Well PCR plates

1. Remove the lock washers (4) from the short screws in the exchangeable thermoblock.
2. Remove the short screws.
3. Connect the adapter plate (2) to the MTP thermoblock (3) using the long screws (1).
4. Secure the screws from below using the lock washers. The lock washers prevent the screws from falling out of the bores if the exchangeable thermoblock is not fixed on the unit.
5. Place the adapter plate on the device with the MTP thermoblock.
6. Tighten the screws.

### 4.1.4 CombiBox

The CombiBox must not be screwed together.

- ▶ Place the CombiBox on the thermsocket.

## 4.2 Inserting tubes

- ▶ Press the tubes lightly into the holders of the thermoblocks.

To maximize temperature control, transfer the micro test tubes in IsoRacks to the thermoblock. The IsoRack reduces the heat exchange of the thermoblock with the environment. If necessary close the IsoRack cover.

### 5 Maintenance

#### 5.1 Cleaning

Clean the exchangeable thermoblocks immediately if sample fluid enters the bore holes or the surface.

1. Clean with a mild soap solution.
2. Rinse thoroughly with distilled water.

#### 5.2 Disinfection / decontamination

---

**Warning! Health hazard from contaminated accessories.**

1. Follow the instructions in the decontamination certificate. It is available in PDF format on our homepage ([www.eppendorf.com/decontamination](http://www.eppendorf.com/decontamination)).
  2. Decontaminate all the parts you want to dispatch.
  3. Include the fully completed decontamination certificate for returned goods in the package.
- 

- ▶ Select a disinfection method complying with the statutory rules and regulations for your area of application. Use e.g. acohol (ethanol, isopropanol) or alcohol-containing disinfectants.

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## 6 Ordering information/Bestellinformationen

Order No. (International)	Order No. (North America)	Description
5361 000.015 5360 000.011 5362 000.019	022670506 022670522 022670549	<b>Exchangeable thermoblock for 24 tubes</b> complete with IsoTherm Rack and IsoTherm cool pack 0°C 24 × 0.5 ml 24 × 1.5 ml 24 × 2.0 ml
5363 000.012	022670565	<b>Exchangeable thermoblock for MTPs and deepwell plates</b> with lid
5363 007.009	022670573	<b>Adapter plate for 96 × 0.2 ml PCR tubes</b> to insert in blocks for MTPs
5363 000.314	022670468	<b>Exchangeable MTP frame</b>
5364 000.016	022670581	<b>Exchangeable thermoblock for 24 micro test tubes</b> 11 - 11.9 mm diam., height 30 - 76 mm
5309 000 333	5309 000 333	<b>Exchangeable thermoblock for 5 ml tubes</b> 8 × 5.0 ml
5366 000.013 5365 000.010	022670531 022670514	<b>Exchangeable thermoblock for Falcon tubes</b> for 8 × 15 ml for 4 × 50 ml
5367 000.017	022670557	<b>Exchangeable thermoblock for 1.5 - 2 ml Cryo tubes</b>
5368 000.010	022670590	<b>Exchangeable thermoblock for slides</b> holds 4 slides, for hybridization experiments
5368 000.100	022670786	<b>Exchangeable thermoblock for slides DC</b> holds 4 DualChip slides, with clamping plate
5352 100.007	022670603	<b>Exchangeable thermoblock CombiBox for any test tubes</b> interior 13.5 × 9.0 × 6.0 cm (W × D × H)
5355 000.011	022670107	<b>Thermomixer comfort</b> without exchangeable thermoblock
5352 000.010	022670204	<b>ThermoStat plus</b> basic device without exchangeable thermoblock 100 - 240 V / 50 - 60 Hz

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Eppendorf North America, Inc. · 102 Motor Parkway · Hauppauge, N.Y. 11788-5178 · USA