


## New spare parts for Thermomixer Comfort 5355 and Compact 5350

Both devices have been modified.


Please note:

The flange of the old motor was lathed (see Fig 1). As a result, the position of the screw heads that hold the motor was deeper. The flange of the new motor has not been lathed (see Fig 2). Instead, the base plate of the new Oscillating Frame System (OFS) has additional holes to make room for the higher standing screw heads. (See Fig 3). However, the „new“ OFS can also be operated with the old motor.

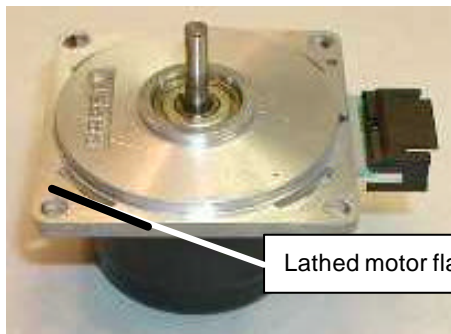


Lathed motor flange

Fig 1 Motor "old version" 5350 005.155

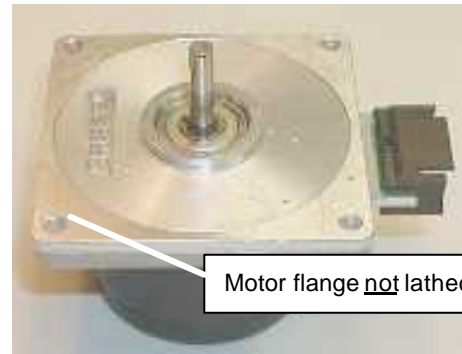

Motor flange not lathed

Fig 2 Motor "new version" 5350 866.002

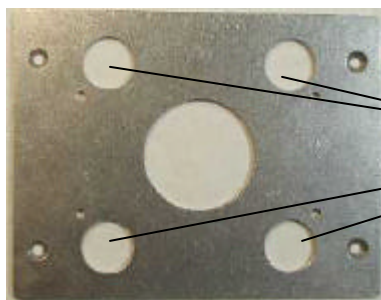

Additional bores for heads of mounting screws of new motor

Fig 3 Base plate OFS "new version".  
View from below.

<b>Motor</b>	<b>5350 866.002 (new)</b>
<b>Oscillating Frame System Comfort</b>	<b>5355 866.000 (new)</b>
<b>Oscillating Frame System Compact</b>	<b>5350 867.009 (new)</b>

Motor 5350 005.155 and OFS 5355 861 009 or 5350 861.000 are still available.

In addition, the shielding in the housing has been changed from red conductive lacquer to silver-colored aluminum damping. The upper part of the housing is compatible and has the order number 5352 853.005.

The heating-cooling plate (HCP) for the Mixer Comfort with brass foil has also been modified. Feature: The sensor of the new HCP is no longer spring-suspended. The HCP is compatible and has the order number 5355 860.002. The order number for the HCP with copper foil 5355 865.004 will be deleted.

The order number for the Thermoblock of the Mixer Compact with brass foil 5350 857.003 will be deleted. The order number for the Thermoblock with copper foil 5350 860.004 has been maintained.

**Chapter 7 of the spare parts list in the service manual has been revised. The manual now has the number 5355 900.055-03.**