

Contact person at HQ, Hamburg:	Ralf Teichmann
Phone:	+494053801-667
E-mail:	teichmann.r@eppendorf.de

New spare part and software version 3.4.3.0 for the BioSpectrometer family 613x and BioPhotometer D30 6133

The spare part “Mirror 6135 866.006” for the BioSpectrometer family “613x” and the BioPhotometer D30 “6133” is no longer available.

The existing mirror will be taken back by Eppendorf until the end of December 2013 and replaced by a new “mirror 6135 866.103“. Please use the SIL procedure for this.



The new mirror will be used in devices from the following serial number onwards:

BioSpectrometer „basic“ Seriennummer: 6135CL601321
 BioSpectrometer „kinetic“ Seriennummer: 6136CJ601387
 BioSpectrometer „fluorescence“ Seriennummer: 6137CJ701337

BioPhotometer D30 Seriennummer: 6133CL400223

If the “mirror 6135 866.006“ is replaced, the new software version 3.4.3.0 must be installed.

In the “Device Information“ service function the article number for the installed mirror must be selected next to “Installed Mirror“. For the new mirror, activate 6135 866.103.

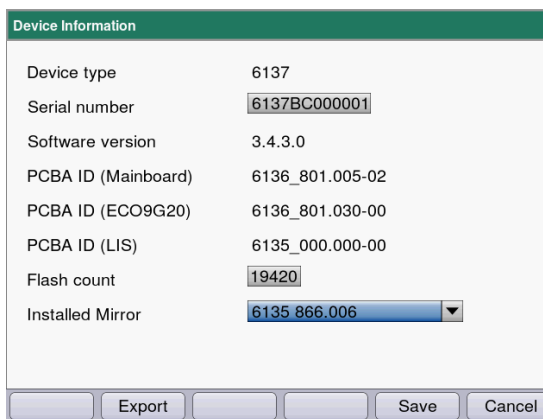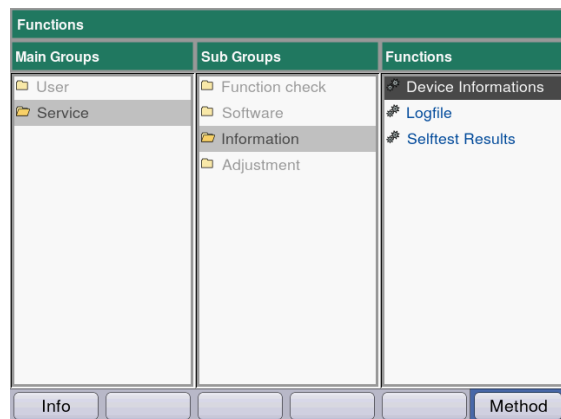

Please note:

- When replacing the main board, these mirror parameters must also be set.
 Existing mirror: activate 6135 866.006
 New mirror: activate 6135 866.103
- After each assembly, the device must be adjusted using the “Optical Adjustment” service function.