

Contact person in HQ, Hamburg:	Ralf Teichmann KS
Phone:	+49 40 53801 667
Email:	teichmann.r@eppendorf.de

## BioSpectrometer (6135, 6136)

### Warranty, transport and repairs

We recommend adding this information to your Service Manual, so that all relevant product information can be found in one place.

We would like to take this opportunity to point out that any warranty claims can only be recognized provided that the product in question has been used in a proper manner in compliance with its intended use. This also applies to the transport conditions for BioSpectrometers.

#### **Always use the original packaging for transport.**

If this packaging has been lost or can no longer be used due to damage, you can request new packaging using the order number 6135 867.002.

BioSpectrometers are sensitive measuring instruments. Is the light path misaligned as the result of improper use (wrong position of the lamp holder), on-site adjustment of the light path will not be possible. It can only be performed at the factory. The Xenon lamp is exchanged by means of a positioning aid (1). This requires that the old lamp holder was correctly positioned.

