

Contact person in HQ, Hamburg:	Ralf Teichmann KS
Phone:	+49 40 53801 667
Email:	teichmann.r@eppendorf.de


## Eppendorf BioSpectrometer® basic / kinetic

A software update (version 1.3.6.0) for the BioSpectrometer is now available as a download at [www.eppendorf-support.com](http://www.eppendorf-support.com).

If the photo diode array (6135 857.007) needs to be replaced during repairs, SW version 1.3.6.0 or a higher version must be installed.

The maximum intensity values in the service function "Intensity Adjustment" have changed with this version. Refer to service manual: **5.5.6.2 Service Function – “Intensity Adjustment”**

- To ensure the correct operation of the BioSpectrometer, peak 1 and the reference sensor must show an intensity of at least 40%.
- The intensity at peak 2 must be a minimum of 29%.

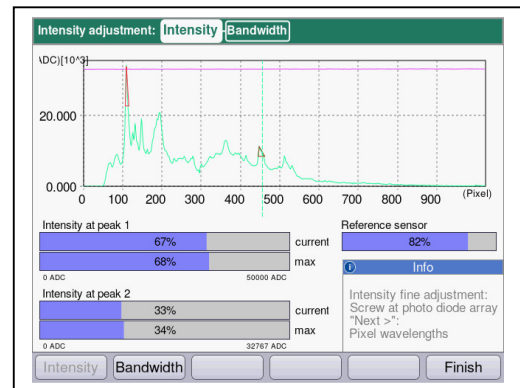

A fault in the help text has also been corrected. The corresponding PACT has been created and will be sent shortly.

We recommend that you add this information to your service manual to ensure that all product-relevant information is filed together in one place.