

Contact person in HQ, Hamburg:	Ralf Teichmann KS
Phone:	+49 40 53801 667
Email:	Teichmann.r@eppendorf.de

BioSpectrometer / BioPhotometer (plus) Certification of Secondary UV-VIS Filter Sets, Test Filter Sets and Reference Filter Sets

We recommend adding this information to your Service Manual.

Product name	<u>Product</u> Order Number International	<u>Product</u> Order Number North America	<u>Certification</u> Order Number International	<u>Certification</u> Order Number North America
BioPhotometer Secondary UV-VIS Filter Set (Customer)	6131 928.007	952010221	6131 928.104	6131928104
BioPhotometer Secondary UV-VIS Filter Set (Service)	6131 931.008	952010239	6131 931.105	6131931105
BioSpectrometer Reference Filter Set (Customer)	6135 928.001	6135928001	6135 928.109	6135928109
BioSpectrometer Test Filter Set (Service)	6135 931.002	6135931002	6135 931.100	6135931100

To ensure correct verification of the photometric and wavelength accuracy of Eppendorf spectrometers and photometers, the filter sets used for this purpose must be checked and recertified every two years. Information on the date on which the last check took place can be found on the certificate which is located on the inside of the filter set box.

Each year EAG offers four fixed dates on which filter sets are checked and certified. This requires that the filter sets concerned are forwarded to EAG via local organizations. Filter sets from end customers also need to be forwarded to EAG with a corresponding order. This order should include the following information: customer order number, certification article number, customer's name and address, customer number, return conditions (air or sea freight, etc.).

Please ensure that all filter sets are dispatched in plenty of time to ensure they reach EAG by the specified dates.

Filter sets which arrive after the date specified can only be checked on the following date!

For 2013 onwards, the following dates apply, unless specified otherwise:

- Start of the second quarter of 2013
- Start of the third quarter of 2013
- Start of the fourth quarter of 2013
- Start of the first quarter of 2014
- and so on....

Delivery address to which the filter sets must be sent:

Eppendorf AG
Goods Receipt – Desinfection
Barkhausenweg 1
22339 Hamburg
Germany

Technical Information



Contact person in HQ, Hamburg:	Ralf Teichmann KS
Phone:	+49 40 53801 667
Email:	Teichmann.r@eppendorf.de

