

Contact person in HQ, Hamburg:	Uwe Krohss
Phone:	+49 40-538 01-193
Email:	Krohss.u@eppendorf.de

Mastercycler ep realplex software under Windows 7/64Bit

The Mastercycler ep realplex software (application software and commissioning software) will not be available as a Windows 7/64 Bit variant.

To ensure that there is a computer capable of running the Mastercycler ep realplex software in the future, the operating system Windows XP/32 Bit or Windows 7/32 Bit is needed.

If it is not possible to provide a computer with the operating system Windows XP/32 Bit or Windows 7/32 Bit, there is the possibility of installing a “virtual machine” (VM) on a computer that is equipped with the operating system Windows 7/64 Bit.

This VM is operated with Windows XP/32 Bit.

It is then possible to install and use all realplex software components inside this “VM” in an unmodified form.

The VMwarePlayer can be downloaded free of charge from

<https://my.vmware.com/web/vmware/details?downloadGroup=PLAYER-312&productId=152>

(after registration).

You also need a Windows XP/32 Bit license.

The following installation guide is also available at www.eppendorf-support.com under “SoftwareDownloads”.