Contact person in HQ, Hamburg:	Uwe Krohss
Phone:	+494053801-193
Email:	Krohss.u@eppendorf.de

## The new nexus Software 2.4.5.0. is available on the Support.com

It is recommended to install this software version to avoid problems with older versions of the software. The new software package also includes the new error list. Should there be any errors after the installation, you can use this error list as a troubleshooting guide, as usual. The first serial number to come out of the EDC with this software version is 6345CO505420 (Lot C2591390).

The software will address and reduce the occurrence of the following errors:

Error #	Description	Change
205	Electrical interferences when setting the temperature	More robust temperature control. The Mastercycler will remain operational even during strong electrical interferences. Check the external power supply in case this error occurs.
408	Error in fan control	Improvement of controller state machine which should prevent this error from occurring.
610	Error in changing state of status LED	Error is ignored. The LED may show a wrong status, e.g. will blink green instead of orange in case of a hold step. This will be fixed with the software for the MC nexus X2.
612	Fan speed to low	The time for initiating the fan control was extended substantially. Should this error still occur, replace the fan first also in together with error 611.
805	Started program could not be found in recently used folder.	This folder shows the last five used programs. Not changed in this software update. This problem will be solved with the software for the MC nexus X2. After closing a program, this program will be removed from that folder.
1412	This message prompts the user to switch the Mastercycler nexus on and off.	This error was introduced to prevent timeout errors 1441 – 1444 from occurring after a fatal error of the temperature control.
1420	The Thermoblock doesn't received an answer from the control unit after starting the program.	The progress bar for the program status is no longer displayed. This will speed up the start-up of the Mastercycler. The error will not occur any more.
1441-1444	Set temperature is not reaching within the corresponding time.	These timeout errors occur as a subsequent error after a fatal error if the Mastercycler has not been switched off and on again. Now it is impossible to start a PCR program after a fatal error has occurred. Error 1412 will now be displayed.
1504	Program is of wrong type for Mastercycler or includes invalid parameters.	The bug which corrupted PCR programs was fixed. This error can only occur if a program is not suitable for the specific Mastercycler model (e.g. tempering rates of a Mastercycler with silver Thermoblock are not executable on a Mastercycler with aluminium Thermoblock).
UK/KS		

Contact person in HQ, Hamburg:	Uwe Krohss
Phone:	+494053801-193
Email:	Krohss.u@eppendorf.de

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