

Contact person in HQ, Hamburg:	Ingo Dietrich
Telefon:	+ 49 40 - 538 01 - 314
Email:	dietrich.i@eppendorf.de

For PCB Power supply unit (6321 856.002) a modified PCB SMPS (6321 895.008) is available.

For Mastercycler pro with PCB Power supply 6321 856.002 the PCB SMPS (switch mode power supply) must be exchanged, if the errors D2C and/or D2D occasionally occur.

For Mastercycler ep 5341 / 5344 / 5345 and Mastercycler realplex 6300 / 6302 the previous PCB Power supply 6321 856.002 as well as the PCB Power supply 5341 856.002 are further on usable without any modifications.


We recommend that you include this information together with Chapter 2 of your Mastercycler Service Manual, to have all relevant product information available in one place.

Für die LP-Netzteil (6321 856.002) ist eine veränderte LP-SMPS (6321 895.008) verfügbar:

Beim Mastercycler Pro mit der LP Netzteil 6321 856.002 muss die LP SMPS (Switch Mode Power Supply) ausgetauscht werden, wenn gelegentlich die Fehler D2C und/oder D2D auftreten.

Im Mastercycler ep 5341 / 5344 / 5345 und im Mastercycler realplex 6300 / 6302 kann sowohl die bisherige LP-Netzteil (6321 856.002) als auch die LP-Netzteil (5341 856.002) weiterhin ohne Änderung eingesetzt werden.

Wir empfehlen, diese Info Ihrem Service Manual Mastercycler Kapitel 2 beizufügen, damit alle relevanten Produktinformationen zusammenhängend abgelegt sind.

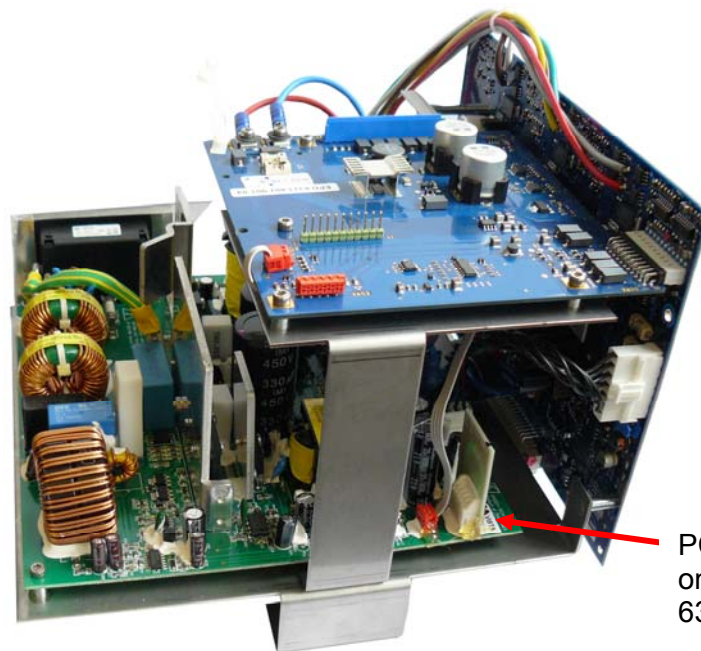

PCB SMPS 6321 895.008
on PCB Power supply
6321 856.002