

We recommend that you include this information together with Chapter 2 of your Service Manual for the Thermomixer/Thermostat, so that all relevant product information can be found in one place.

Thermomixer compact 5350
Thermomixer comfort 5355
Thermostat plus 5352

Adapter for temperature adjustment
Adapter cable for temperature adjustment 5350

The adapter for temperature adjustment and the adapter cable for temperature adjustment 5350 for automatically adjusting the temperature of the three devices can no longer be produced. These two auxiliary aids are no longer available. The temperature adjustment can now only be performed manually. A description of how to adjust the temperature manually can be found in the Service Manual (Chapter 7).

In order to facilitate the temperature measurement for manual temperature adjustment, and to minimize measurement errors, we have developed a temperature sensor in the shape of a 1.5 mL tube. Use this temperature sensor in combination with the temperature measuring device from the temperature validation system MC / MC ep (0055 000.298) for manually adjusting the temperature on all three devices.

The temperature measuring device and the temperature sensor 1.5 mL can also be used for temperature calibration of the devices. Further information on this topic can be found in the operating manual for the temperature sensor 1.5 mL and for the temperature validation system MC / MC ep.

Order numbers **old** (no longer available):

0024 204.707 Adapter for temperature adjustment

0024 204.804 Adapter cable for temperature adjustment 5350

Order numbers **new**:

5354 850.500 Temperature sensor 1.5 mL

(0055 000.298 Temperature validation system MC / MC ep)